

# Microstructural Investigation of Austempered Ductile Iron (ADI) with “Shuttle & Find” Interface for Correlative Microscopy in Materials Analysis

Authors: C. Scherrer, Institute of Materials and Process Engineering, ZHAW, Winterthur, Switzerland  
D. Lysenkov, H. Mantz, Carl Zeiss NTS GmbH, Oberkochen, Germany  
C. Thomas, Carl Zeiss AG, Corporate Research & Technology, Oberkochen, Germany

Date: November 2009

## Introduction

Austempered Ductile Iron (ADI) excels through strength, wear resistance and toughness – characteristics that make ADI the material of choice for use in combustion engines and gear box components. This means that safety aspects are also involved in addition to purely functional aspects. For this reason, changes in the ADI production process need to be monitored with respect to the material's characteristics and must be optimized systematically. For the micro- and nanoscopic analysis of the structure and precipitations, scientists typically use both light and electron microscopes. To date, however, there has been no possibility of relocating regions of interest without doubt when transferring the sample from the light to the electron microscope or vice versa. “Shuttle & Find” – the interface for correlative microscopy in materials analysis offers an easy-to-use solution, allowing seamless integration of these two complementary technologies for the first time.

## Material

EN-GJS-1200-2 according to DIN EN 1564.

Austempered Ductile Iron (ADI) is a family of iron-based materials, which among different processes can be obtained by heat treatment (so-called austempering) of nodular cast iron. ADI is well-known for its excellent strength, wear resistance and toughness. The ultimate tensile strength of min. 1200 MPa and elongation at fracture of min. 2 % results in a cost-effective solution providing comparable performance to high strength aluminum alloys (per unit weight) or even steels.




## Equipment

The current investigations were done on a ZEISS Axio Imager.M1 light microscope with motorized stage, SUPRA™ 40VP Field Emission-SEM (FE-SEM) equipped with an AsB® detector, SUPRA™ 55VP FE-SEM with an AsB® and Bruker Quantax 200 EDS detectors.

In general, in order to enable the correlative workflow, the following points should be fulfilled:

- Light microscope equipped with digital camera and motorized stage of Axio Imager, Axio Observer and SteREO Discovery families with corresponding mounting frames,
- Scanning electron microscope of EVO®, SIGMA, SUPRA™, ULTRA and MERLIN® families with an SEM adapter for correlative microscopy,
- Correlative holder,
- Software compatibility requirements.

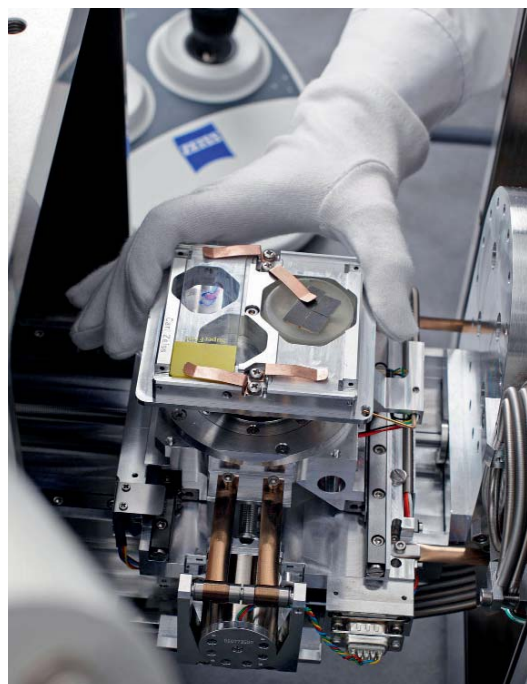

## Application

Supporting the tribological characterization the main investigation task was an exact description, including elemental analysis, of the materials microstructure. This was, however, only possible with the help of correlative microscopy, because the (hard) precipitates allowed a systematic investigation in the scanning electron microscope (SEM) after being examined in the light microscope only in case of a precise relocation. In addition, the task became more complicated by the fact that it was only possible to detect the precipitates in the SEM with the backscatter electron detector (e.g. AsB®).

The identification of the chemical composition of precipitates is possible only by EDS. Thus, the exact relocation of the region of interest (ROI) – in order to avoid time consuming searches – is of great importance. “Shuttle & Find” – correlative microscopy solution from Carl Zeiss fulfills exactly these application requirements and allows micron-precise relocation of the ROI on metallographic samples even at high magnifications in both light and electron microscopes, making subsequent EDS in a SEM a matter of routine (Figs. 1–4).

## Results

The matrix basically consists of bainite (in fact ausferrite), graphite nodules and of individual areas of retained austenite. The precipitates were found in these areas. After the relocation of the precipitates in SEM with the help of “Shuttle & Find”, the area was scanned with EDS for Mo, Fe, P and C. The spatial resolution of EDS in this experiment is limited in order to provide information about the exact position of the neighboring elements. Nevertheless, the EDS mappings showed that Mo, P and C are found in the same area, verifying the existence of iron phosphides and molybdenum carbides respectively. The wear resistance increases through such hard precipitates whereas the toughness decreases.

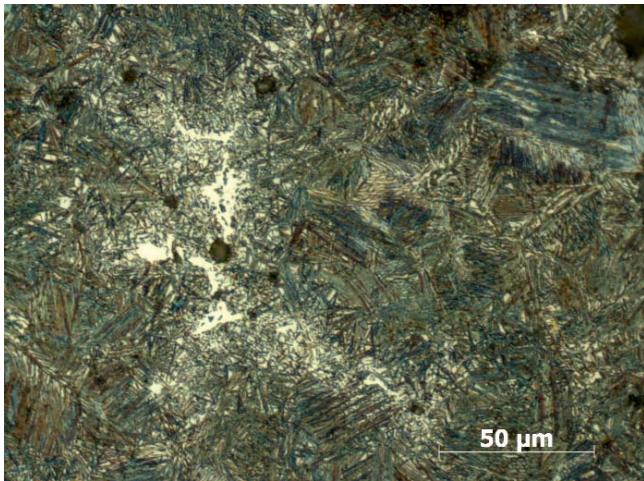


Fig. 1. Light microscope image of ADI sample with a magnification of approx. 400:1. One can see the ROI with the precipitates.

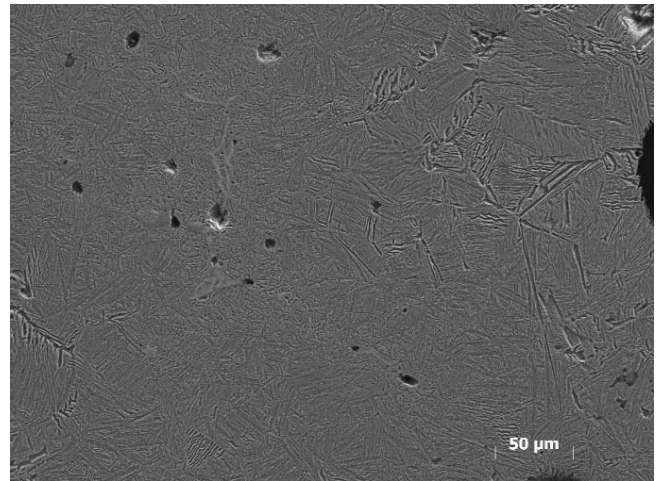


Fig. 2. SE image of the same area as in Fig. 1. The problem of relocating the same area with SE becomes evident.

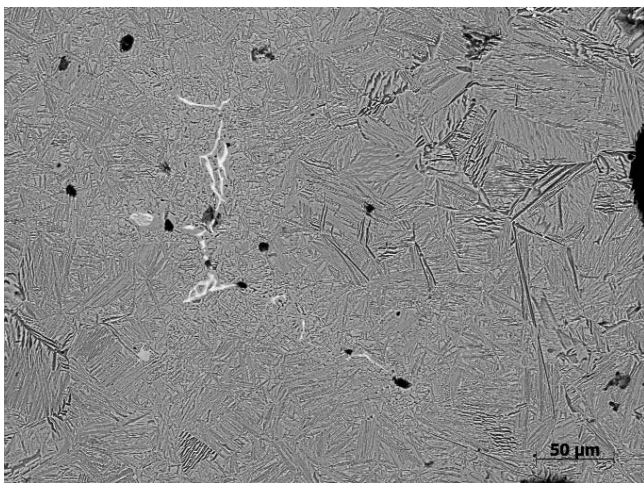


Fig. 3. BSE image of the same area as in Fig. 1. The microstructure is clearly visible in comparison to the SE image.

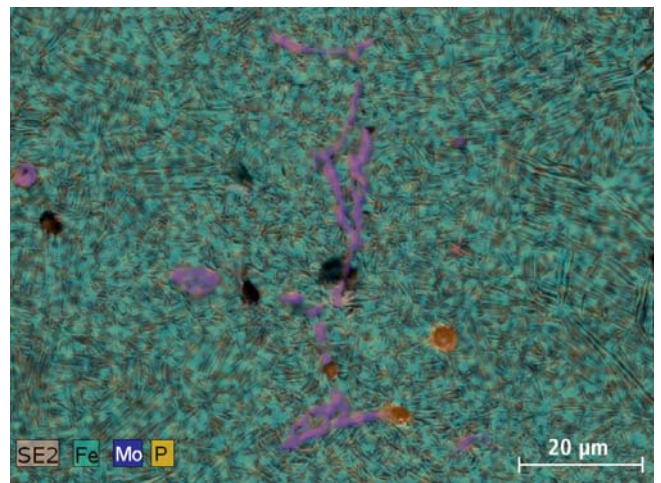


Fig. 4. SE image of the same area as in Fig. 1 overlapped with Mo, Fe and P mappings obtained with EDS.

Note: The Mo (La) line (2.29 keV) and S (Ka) line (2.31 keV) interfere. An acceleration voltage of 30 kV and a long analysis time are required in order to resolve Mo.

**Acknowledgements:**

We thank the Institute of Materials and Process Engineering, ZHAW, Winterthur (Switzerland) for providing the samples.

**Maximum Information – Maximum Insight**

More than 160 years of experience in optics has laid the foundation for pioneering electron and ion beam microscopes from Carl Zeiss. Superior integration of imaging and analytical capabilities provides information beyond resolution, unlocking the best kept secrets of your sample.

With a broad technology portfolio Carl Zeiss provides instruments both tailored to your requirements and adaptable to your evolving needs. With our highly versatile application solutions we endeavor to be your partner of choice.

Superbly equipped, regional demo centers provide you with access to our applications expertise developed in collaboration with world-class partners in industry and academia. Global customer support is provided by the Carl Zeiss group together with an extensive network of authorized dealers.

Our mission at all times: Maximum Information – Maximum Insight.

**Carl Zeiss NTS GmbH**

A Carl Zeiss SMT AG Company  
Carl-Zeiss-Str. 56  
73447 Oberkochen, Germany  
Tel. +49 73 64 / 20 44 88  
Fax +49 73 64 / 20 43 43  
info-nts@smt.zeiss.com

**Carl Zeiss SMT Inc.**

One Corporation Way  
Peabody, MA 01960, USA  
Tel. +1 978 / 826 1500  
Fax +1 978 / 532 5696  
info-usa@smt.zeiss.com

**Carl Zeiss SMT Pte Ltd.**

50 Kaki Bukit Place #04-01  
Singapore 415926, Singapore  
Tel. +65 65 67 / 30 11  
Fax +65 65 67 / 51 31  
info.sea@smt.zeiss.com

**Carl Zeiss MicroImaging GmbH**

Königsallee 9-21  
37081 Göttingen  
Germany  
Tel. +49 551 50 60 660  
Fax +49 551 50 60 464  
micro@zeiss.de  
[www.zeiss.de/micro](http://www.zeiss.de/micro)  
[www.zeiss.de/correlative\\_microscopy](http://www.zeiss.de/correlative_microscopy)

**Carl Zeiss SMT Ltd.**

511 Coldhams Lane  
Cambridge CB1 3JS, UK  
Tel. +44 12 23 41 41 66  
Fax +44 12 23 41 27 76  
info-uk@smt.zeiss.com

**Carl Zeiss SMT S.a.s.**

Zone d'Activité des Peupliers  
27, rue des Peupliers -  
Bâtiment A  
92000 Nanterre, France  
Tel. +33 1 41 39 92 10  
Fax +33 1 41 39 92 29  
info-fr@smt.zeiss.com

[www.smt.zeiss.com/nts](http://www.smt.zeiss.com/nts)

[www.smt.zeiss.com/correlative-microscopy](http://www.smt.zeiss.com/correlative-microscopy)



We make it visible.